Version : 2024 Rev.0

VACUUM PUMP

V Series

SH-V10/V20/V40/V60/V100

TABLE OF CONTENTS

Thank you for purchasing our product. We know that in today's competitive marketplace, customers have many choices when purchasing laboratory equipment. We appreciate your choosing our quality product. We stand behind our products and want to let you know we are here if you need us. Before you use the unit, read this entire manual carefully to understand how to install, operate and maintain the unit in a safe manner. Your satisfaction with the unit will be maximized as you read this manual thoroughly. Our capable products will satisfy you by the best performance with easy operation.

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| | |
|-----------------------------------|----|
| 1. Features & Specification | 3 |
| 2. Basic Parts | 5 |
| 3. Installation | 6 |
| 4. Operation | 7 |
| 5. Inspection & Maintenance | 8 |
| 6. Troubleshooting | 12 |
| 7. Warranty | 15 |

1. Features & Specification

SH-V Series

Perfect design and performance

Constantly pursuing on highest technology norms, SH Scientific V series vacuum pump reached new highest level: cylinder structure and innovative oil circulation system results in super stability; auto anti-suck back valve and two shifts gas ballast makes best reliability; Perfect combination of technology and aesthetics helps to become classic.

- Tow-Shift adjustable gas ballast valve satisfies different requirements of condensable vapor (such as water vapor) to be exhausted out of pump in different processes.
- Dual protection of oil anti-suck back ensures vacuum system from oil pollution when pump stops running and needs to be easily restarted.
- Forced oil circulation system consisted of oil pump and constant pressure oil supply mechanism ensures stable running of the pump.
- Less components are used, easy to maintain and repair.

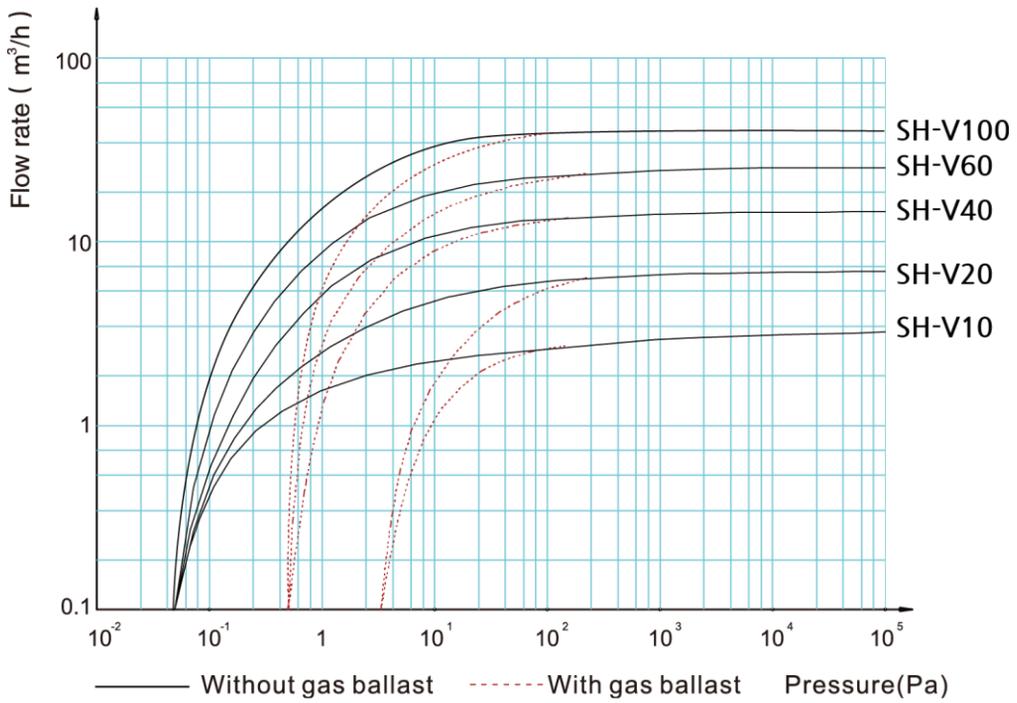
Specifications

| Model | SH-V10 | SH-V20 | SH-V40 | SH-V60 | SH-V100 |
|---|--|--|--|--|--|
| Pumping Speed(60Hz) | 78L/min | 156L/min | 312L/min | 474L/min | 960L/min |
| Ultimate total pressure gas ballast close | 5x10 ⁻¹ Pa (3.7x10 ⁻³ torr) | 5x10 ⁻¹ Pa (3.7x10 ⁻³ torr) | 4x10 ⁻¹ Pa (3.0x10 ⁻³ torr) | 4x10 ⁻¹ Pa (3.0x10 ⁻³ torr) | 4x10 ⁻¹ Pa (3.0x10 ⁻³ torr) |
| Ultimate total pressure gas ballast open | 10 Pa (7.5x10 ⁻² torr) | 10 Pa (7.5x10 ⁻² torr) | 8x10 ⁻¹ Pa (6x10 ⁻³ torr) | 8x10 ⁻¹ Pa (6x10 ⁻³ torr) | 8x10 ⁻¹ Pa (6x10 ⁻³ torr) |
| Power supply | 220V, 1Φ, 50/60Hz | | | | 220V, 3Φ, 50/60Hz or 380V, 3Φ, 50/60Hz |
| Power rating(kW) | 0.4 | 0.4 | 0.75 | 1.1 | 1.5 |
| Inlet/Exhaust Flange | KF 16/25 | KF 16/25 | KF 25 | KF 25/40 | KF 40 |
| Oil capacity(L) | 0.6~1.0 | 0.6~1.0 | 0.9~1.5 | 1.3~2.0 | 3.3~4.5 |
| Ambient temperature(°C) | 10~40°C | | | | |
| Noise level(dB) | ≤56 | ≤56 | ≤58 | ≤58 | ≤62 |
| Weight(kg) | 19 | 19 | 30 | 35 | 62 |


Option: Oil Mist Trap

Flow Rate Characteristic


Dimension

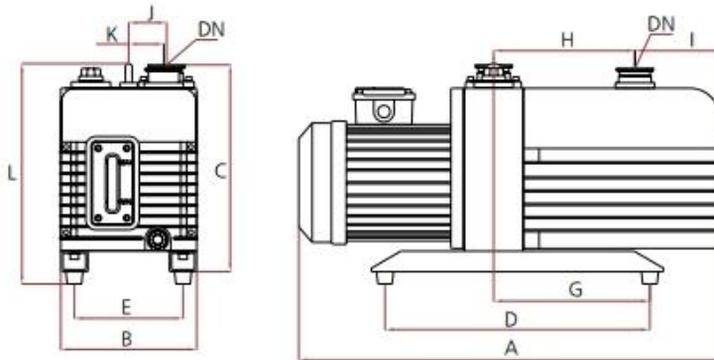Dimensions


| Model | A | B | C | D | E | G | H | I | J | K | L | DN |
|---------|-----|-----|-----|-----|-----|-----|-------|-----|----|----|-----|---------|
| SH-V10 | 440 | 144 | 217 | 210 | 110 | 95 | 143.5 | 45 | 45 | 34 | / | KF16/25 |
| SH-V20 | 440 | 144 | 217 | 210 | 110 | 95 | 143.5 | 45 | 45 | 34 | / | KF16/25 |
| SH-V40 | 530 | 188 | 272 | 320 | 148 | 160 | 165 | 69 | 59 | 38 | 295 | KF25 |
| SH-V60 | 567 | 188 | 272 | 320 | 148 | 160 | 165 | 69 | 59 | 38 | 295 | KF25 |
| SH-V100 | 730 | 234 | 358 | 396 | 190 | 200 | 223 | 157 | 69 | 55 | 390 | KF40 |

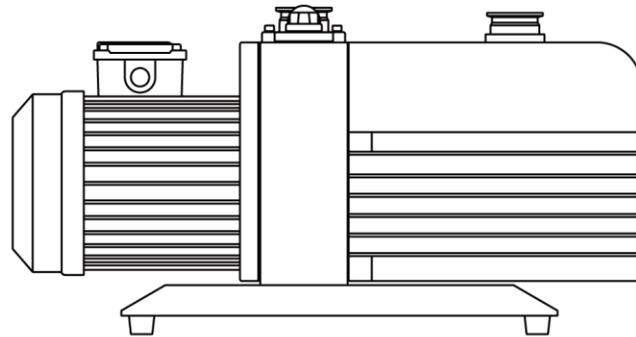2. Basic Parts


<Main unit>

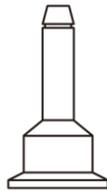+


Hose barb : 1EA

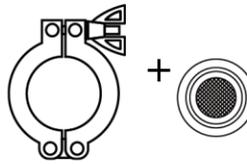Clamp with O-ring: 2SETS

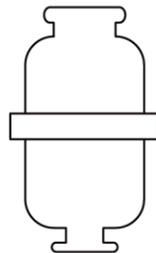
Oil


Oil Mist Trap(Optional)

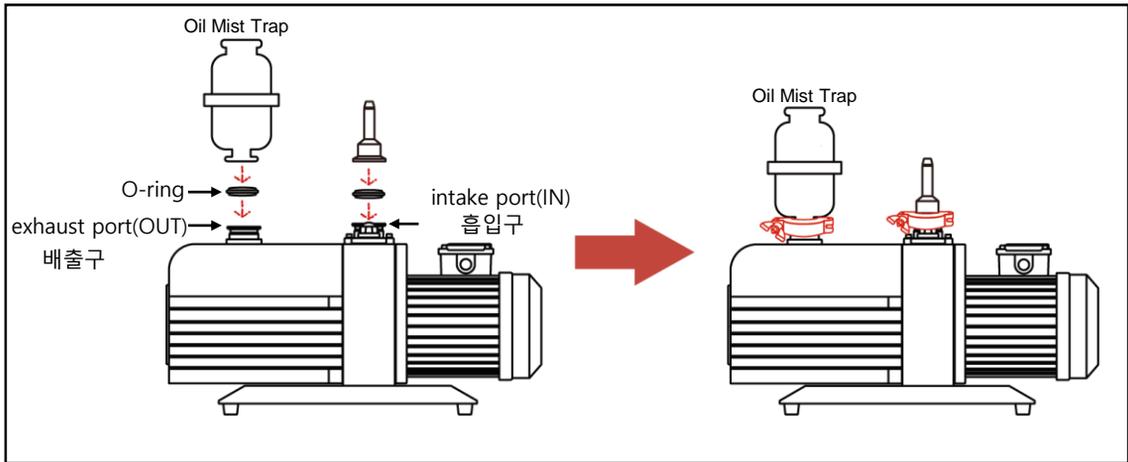3. Installation

- ① Please put the O-ring between the intake port and the hose barb and tighten the clamp.
 (*In case of Oil Mist Trap, please install it in the exhaust port in same way.)

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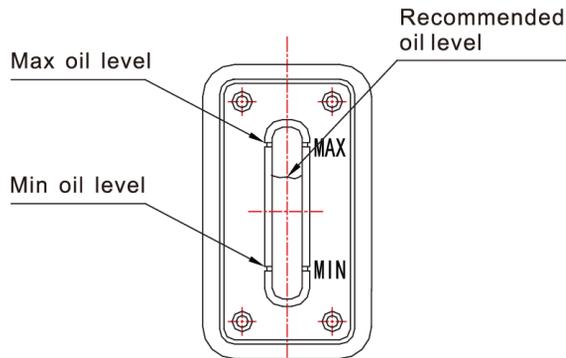
<Fig. 1>

- ② The oil level shall be kept between position MAX and MIN of oil sight glass.
 Excessive or insufficient oil will decrease the pump performance or even cause malfunction. The pump must be stopped before adding oil.

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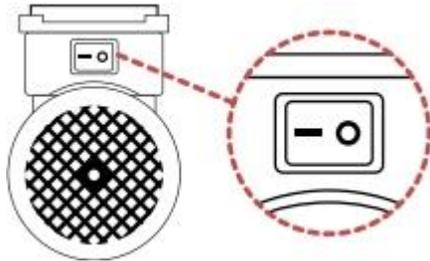4. Operation

- ① Please plug-in the power cord 220V and switch on the pump.

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Gas ballast valve knob arrow

<Fig. 2> Gas Ballast Valve

- ② Please lock the gas ballast valve. (Set to the position C, Fig.2)

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High temperature hazard! Do not touch the surface! The surface of the pump could be very hot during operation and after shutdown within one hour.

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Make sure the gas flow at the exhaust port is not blocked or restricted in anyway. Do not start the pump if the exhaust port is blocked.

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The operating ambient temperature is 10~40°C and humidity should be less than 85%.

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5. Inspection and Maintenance

All checks and maintenance must be performed by trained personnel in accordance with local safety rules and regulations.

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In case of presence of hazardous substances due to oil decomposition or media sucked into the pump, it's necessary to define the hazard and take all necessary safety precautions.

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
In the presence of hazardous substances, determine the hazard in the first time and follow all appropriate safety procedures. If the potential hazard still exists, the pump must be decontaminated before any maintenance.

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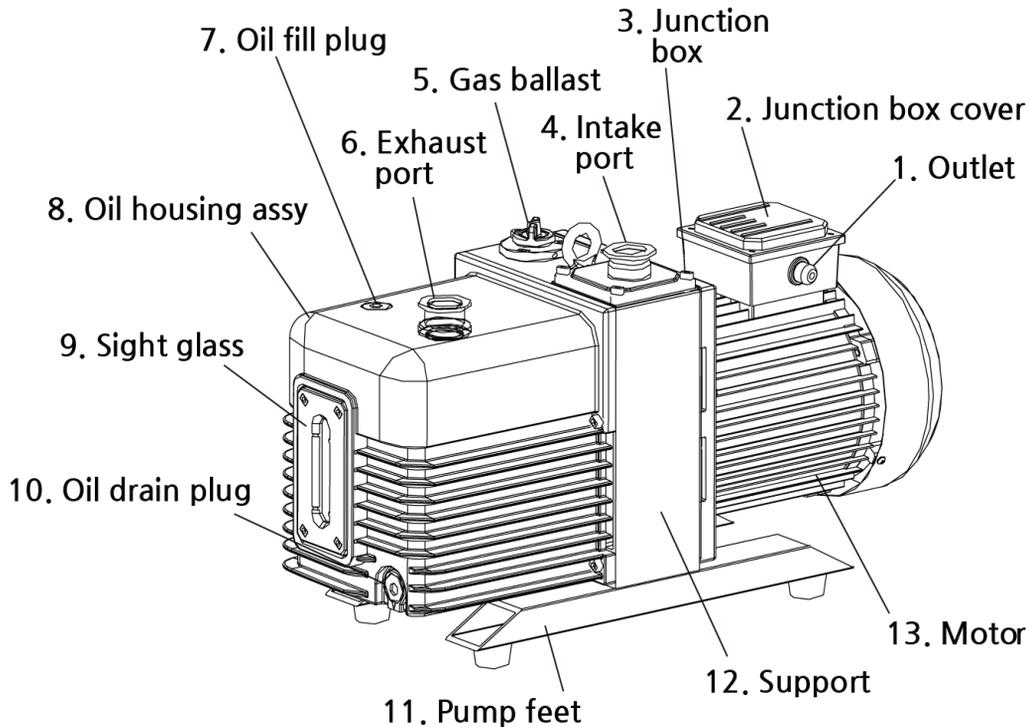<Fig. 3> Outside view

• 5.2.1 Checking oil level 오일레벨 확인

- 1) The oil level shall be always kept between position MAX and MIN of oil sight glass during operation (refer to Fig. 1) Add oil when the oil level is lower than position MIN and discharge oil by removing the drain plug (refer to Fig. 3) when the oil level is above position MAX.

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• 5.2.2 Checking pump's noise 펌프소음 확인

The pump noise shall be continuous and stable during operation. If any abnormal noise is found, please check troubleshooting.

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- 1) Change the oil when the pump is switched off and cooled down.
전원을 끄고 펌프를 식힌 후, 오일교체하십시오.
- 2) Remove the oil drain plug(Fig.3) and drain the used oil into a suitable receptacle. When the flow of oil stops, screw on the oil drain plug and start the pump shortly(10 seconds). Switch off the pump, remove the plug again and drain the residual oil. Screw on the oil drain plug (Check O-ring, replace it if necessary). Remove the oil fill plug(Fig.3) and fill in fresh pump oil which is specified or provided by the manufacturer.
(Fig. 3) 오일배유 플러그를 제거하여 사용한 오일을 배출시킵니다. 배출되는 오일 흐름이 멈추면, 오일 배유플러그를 잠그고 10초 정도만 펌프 작동 시킵니다. 그런 후, 펌프 전원을 끄고 오일 배유 플러그를 다시 제거하여 잔여 오일을 배출시킵니다. 다시 오일 배유플러그를 잠그고(O-ring을 체크하고 필요시 교체), (Fig. 3)처럼 오일보충플러그를 제거하여 펌프제조사 지정된 새로운 오일로 교체하십시오.
- 3) After oil changing, handle the used oil stored in an assigned container according to local relevant environmental laws and regulations.
오일 교체 후, 사용한 오일은 안전수칙에 맞게 적합한 용기에 보관 관리 하십시오.

• 5.2.4 Regularly checking intake filter and anti-suck back valve

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Regularly take down the intake prot and clean the filter by compressed air. Check the seal position of anti-suck back valve for cleanness, damage, hardening and height.

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• 5.2.5 Regularly checking fan cover and motor fan

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Regularly remove the clean the fan cover and motor fan. Clean them by compressed air and reinstall them properly before pump operation.

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7. Warranty

SH-V series vacuum pump have one year guarantee from the date of purchase. Spare parts will be provided free of charge within the period of guarantee in condition that the pump is operated according to the operating manual.

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The following failures will be charged for repair:

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E-mail. sales@samheung21.com

1544-0351

Overseas sales office www.samheung21.com

Daejeon, Republic of Korea

Tel. 042) 826 5172 Fax. 042) 826 5173

E-mail. i-sales@samheung21.com

SH Scientific Corporation USA labandfurnace.com

Portland, Oregon

E-mail. jhbang@samheung21.com