

Version : 2021 Rev.0

Dual-Stage Low noise rotary
VACUUM PUMP



VDC (Low noise) Series
SH-VDC7 / VDC10 / VDC20

TABLE OF CONTENTS

Thank you for purchasing our product. We know that in today's competitive marketplace, customers have many choices when purchasing laboratory equipment. We appreciate your choosing our quality product. We stand behind our products and want to let you know we are here if you need us. Before you use the unit, read this entire manual carefully to understand how to install, operate and maintain the unit in a safe manner. Your satisfaction with the unit will be maximized as you read this manual thoroughly. Our capable products will satisfy you by the best performance with easy operation.

1. Use Information	3
2. Attention	4
3. Description	6
4. Technical Parameters	7
5. Transport and Storage	8
6. Installation and Connection	8
7. Initial Start and Operation.....	10
8. Inspection and Maintenance	11
9. Troubleshooting.....	13
10. Warranty.....	15

1. Use Information

Dear distinguished customer,

Thanks for choosing our vacuum pump of high reliability (Hereinafter referred to as "the pump").

Please check carefully whether the product received is the same as you ordered and the accessories, spare parts & operating manual are attached as well. Please also check if there's any damage occurred during transportation. If needed, contact the local distributor or our sales team.

In order to maintain a stable performance level of the pump, read this operating manual carefully to fully understand the safety instructions, technical data as well as operating procedures before installation, operation, repair and maintenance of the pump.



Warning

Failure to observe the terms could result in serious personal injury.



Notice

Failure to observe the terms could result in damage to the pump.



This warning label indicates risk of electrical shock. Disconnect the pump from the power supply before beginning with connections, repair and maintenance. Make sure the cover of junction box is properly installed before running.



This warning label indicates high temperature hazard. Do not touch the pump when the pump is in operation.



Notice

Read the operating manual carefully and follow the operating procedures. We reserve the right to modify the design and technical data of the pump without notice which may have discrepancies in the manual. Add vacuum oil as requested before starting the new pump.

2. Attention

In order to ensure the personal safety, read the operating manual carefully before installation, operation, repair and maintenance.



Warning

According to the technical guidelines and wiring regulations for electrical equipment, the power supply must match with the information shown on the pump nameplate and the electrical connections must be performed by a skilled electrician.



Warning

To prevent electrical hazards and sudden start-up hazard which may result in personal injury, disconnect the power supply before checking or repairing the pump.



Warning

Before starting the pump, the motor must be effectively grounded and properly connected with a rated motor protection switch.



Warning

The pump is strictly prohibited to pump dusty, active toxic, corrosive, flammable and explosive gases.



Warning

Do not place obstacles which may affect the ventilation of the motor to avoid burn hazard and fire caused by abnormal temperature rise.



Warning

The exhaust passage must be unimpeded during operation. Make sure that the gas flow from the exhaust port is not blocked or restricted in any way.



Notice

The pump must be operated at ambient temperature between 10-40°C



Notice

Check the oil level before running. Do not operate the pump without oil or short of oil. Otherwise it will result in pump failure.



Notice

Check oil leakage of the pump to prevent slip hazard caused by spilled oil.



Notice

Running without oil lubrication or wrong rotation direction of the motor may cause pump failure.

**Notice**

Do not connect the intake port with an overpressure device. Make sure the size of the exhaust passage with absolute pressure is not more than 1.35 bar (relative pressure is not more than 0.35 bar).

**Notice**

Disposal of used oil and other parts must be in compliance with local relevant environmental laws and regulations.

**Notice**

To pump condensable gases, only use the pump installed with gas ballast.

**Notice**

While pumping a small amount of dust and condensable gases, an installation of corresponding filters is required. Otherwise it will cause pump failure or sharp drop of performance.

3. Description

VDC series vacuum pump is an innovative motor direct drive dual-stage oil sealed rotary vane pump with DC motor. The pump features small motor power design for energy-saving, DC output with low noise level and optional corrosion resistance. Therefore, the pump is an ideal choice for high requirements of quiet operation, especially for analytical applications.

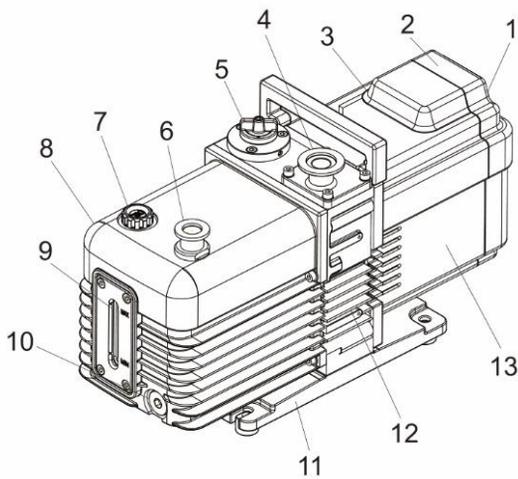


Fig. 1 Outside view

1. Outlet
2. Junction box cover
3. Junction box
4. Intake port
5. Gas ballast
6. Exhaust port
7. Oil fill plug
8. Oil housing assy
9. Sight glass
10. Oil drain plug
11. Pump feet
12. Support
13. Motor

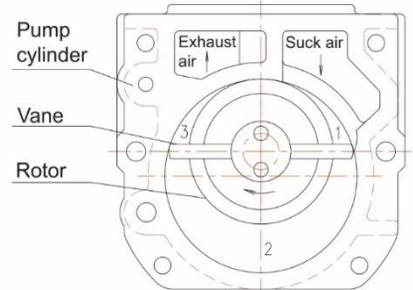


Fig. 2 Functional diagram

3.1 Principle of Operation

Refer to Fig. 2, the rotor mounted eccentrically in the pump cylinder, has two vanes which divide the pump chamber into three changeable compartments. When the rotor rotates clockwise, gas is sucked into chamber 1 gradually, transferred by chamber 2, and compressed and exhausted by chamber 3. The working cycle is repeated to achieve the vacuum of the system. The exhausted gas from the first stage goes into the second stage (low-vacuum stage) through air channel, and is compressed again before discharging into atmosphere to achieve a lower ultimate vacuum pressure.

3.2 Principle of Gas Ballast

The controllable gas (generally dry air at room temperature) is admitted into the pump chamber. It is mixed with pumped vapor during the gas compression process. This increases the percentage of non-condensable gas such that the partial pressure of the vapor being pumped is below its saturated vapor pressure when the exhaust valve opens. The vapor is then discharged from the pump without liquefaction. The more vapor contained in the pumped gas, the more dry air is required (See Fig. 7 Gas ballast valve).

4. Technical Parameters

Model	SH-VDC7	SH-VDC10	SH-VDC20
Pumping Speed(50/60Hz)	66L/min	96L/min	183L/min
Ultimate total pressure gas ballast close	5×10^{-2} Pa (3.7×10^{-4} torr)	5×10^{-2} Pa (3.7×10^{-4} torr)	5×10^{-2} Pa (3.7×10^{-4} torr)
Ultimate total pressure gas ballast open	100 Pa	100 Pa	100 Pa
Power supply	220V,50/60Hz or 110V,50/60Hz		
Power rating(kw)	0.37(1 Phase)	0.37(1 Phase)	0.37(1 Phase)
Inlet/Exhaust Flange	KF 16/25	KF 16/25	KF 16/25
Oil capacity(L)	0.6~1	0.6~1	0.6~1
Ambient temperature(°C)	5~40		
Noise level(dB)	≤50	≤50	≤54
Weight(kg)	14.5	14.5	15.5

Table 1

Pumping speed characteristics

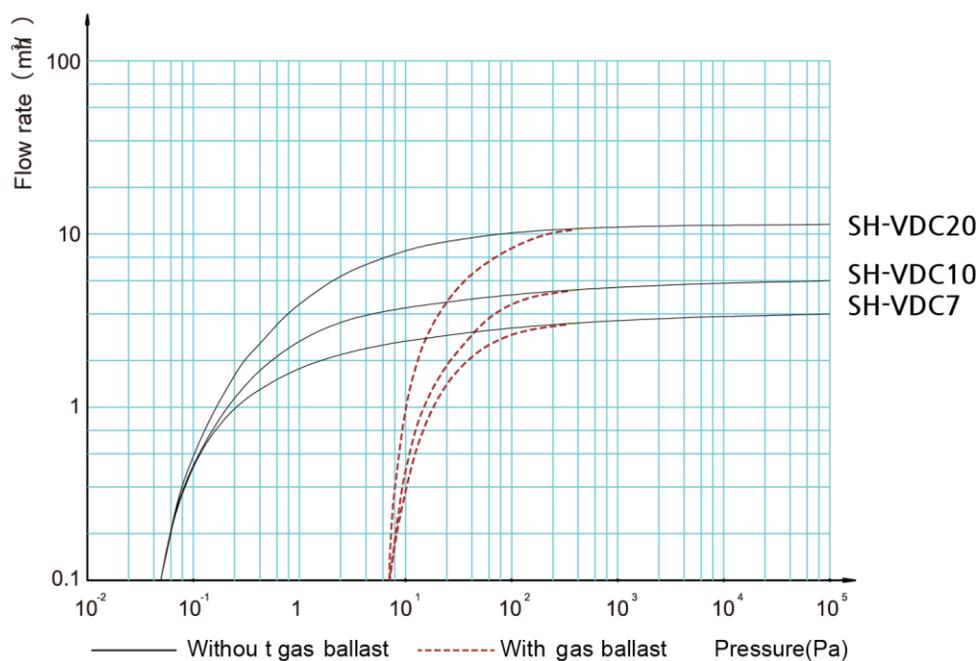


Fig. 3

5. Transport and Storage



Warning

Do not move the pump unless it's switched off and the power supply is disconnected.



Notice

Check carefully if there's any packaging damage during transportation.



Notice

Move the pump with care. Any negligence will cause pump damage during transportation. In order to avoid oil spilling, move the pump fully filled with oil in horizontal direction.



Notice

Disposal of used packing material shall be in compliance with local environmental laws and regulations.



Notice

For long term storage, clean the pump and empty the oil inside. Store the pump in a dry place at room temperature.

6 Installation and Connection

6.1 Installation Dimensions

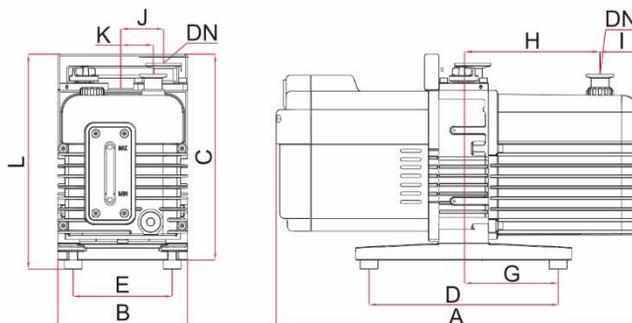


Fig. 4

Dimensions												
Model	A	B	C	D	E	G	H	I	J	K	L	DN
SH-VDC7	420	144	207	210	110	126	154	45	45	34	/	KF16/25
SH-VDC10	420	144	207	210	110	126	154	45	45	34	/	KF16/25
SH-VDC20	420	144	207	210	110	126	154	45	45	38	/	KF16/25

Table 2

6.2 Connection to System



Warning

The pump is strictly forbidden to operate in areas with risk of fire and explosion.



Warning

Make sure the gas flow at the exhaust port is not blocked or restricted in any way. Check to ensure the exhaust line is not obstructed by deposits before start-up.



Warning

The exhaust pressure shall not be 1.15 bar higher than the absolute pressure (relative pressure 0.15 bar) but not be lower than atmospheric pressure.



Notice

Take the following into account to choose the pump installation site: convenient for installation, maintenance, disassembly and electrical connections as well as good ventilation.



Notice

Place the pump feet horizontally or fix the feet with bolts while connecting the pump to vacuum system.



Notice

Oblique installation may result in pump's vibration, high noise or even damage. The pump shall be set up on a flat and firm surface.



Notice

The length of connecting line between pump and vacuum system shall be as short as possible. The diameter of the connecting line shall be the same as that of intake port.



Notice

The hoop at intake port is available for connection with vacuum hose or regular vacuum pipe. No any stress generated by the pipe on the hoop is allowed while connecting.



Notice

The exhaust line should be laid so that it slopes down so as to prevent condensate from backstreaming into the pump and contaminating the oil. Drain the oil in the exhaust line periodically to avoid block of exhaust line. If the exhaust line slopes up, the installation of a condensate trap is necessary. Check leak of the connection between pipe and flange. Vacuum tightness is essential for the achievement of ultimate pressure of the pump.

6.3 Power Supply



Warning

Make sure the power supply is correct. The electrical connections must be performed by a skilled electrician according to the technical guidelines and wiring regulations for electrical equipment. The extra selected cable and motor protection switch shall have the same rated current value as stated on the motor nameplate.

7. Initial Start and Operation

7.1 Start-up



Warning

Make sure the gas flow at the exhaust port is not blocked or restricted in any way. Do not start the pump if the exhaust port is blocked.



Warning

The oil level shall be kept between position MAX and MIN of oil sight glass. Excessive or insufficient oil will decrease the pump performance or even cause malfunction. The pump must be stopped before adding oil.



Warning

Make sure the gas flow at the exhaust port is not blocked or restricted in anyway. Do not start the pump if the exhaust port is blocked.

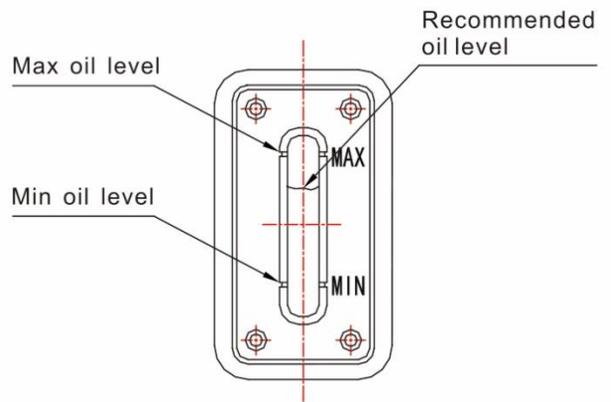


Fig. 5



Notice

The operating ambient temperature is 10 - 40°C and humidity should be less than 85%.

7.2 Operation



Notice

High temperature hazard! Do not touch the surface! The surface of the pump could be very hot during operation and after shutdown within one hour.

Pumping non-condensable gases

Close the gas ballast to pump non-condensable gases (Set the gas ballast to position C, refer to Fig. 6). The ultimate pressure will rise (loss of vacuum) with gas ballast opened.

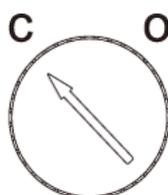


Fig. 6 Gas Ballast Valve

Pumping condensable gases and vapor

If the vacuum system contains a small amount of condensable gas, open the gas ballast valve (Set the gas ballast to position I or II, refer to Fig. 6) and keep the pump running for at least 30 minutes to pump the condensable gas effectively. Close the gas ballast valve when the vacuum system pressure is reduced to a certain value. If the pump operates at low temperature, condensable gas may be dissolved in the pump oil. The oil may be emulsified or deteriorated which will affect the pump performance, even cause corrosion of the pump body.

The rise of oil level during pump operation is a sign of condensable gas contained in the vacuum system. Do not switch off the pump immediately when the process is finished. Keep the pump running with gas ballast open and intake port closed until all gases dissolved in the oil are removed.

8. Inspection and Maintenance



Warning

All checks and maintenance must be performed by trained personnel in accordance with local safety rules and regulations.



Warning

In case of presence of hazardous substances due to oil decomposition or media sucked into the pump, it's necessary to define the hazard and take all necessary safety precautions.



Warning

In the presence of hazardous substances, determine the hazard in the first time and follow all appropriate safety procedures. If the potential hazard still exists, the pump must be decontaminated before any maintenance.

8.1 Routine inspection

Content	Operation	Interval	Remarks
Oil level	Visual inspection	Every three days	5.2.1
Oil color	Visual inspection	Every three days	5.2.1
Pump noise	Acoustic inspection	Every three days	5.2.2
Pump vibration	Visual inspection	Every three days	Check the pump feet or feet bolts in case of abnormal vibration.
Pump temperature	Temperature meter	Every one week	Check and clean fan of the pump and motor for deposits.
Seal & O-ring	Visual inspection	Every one month	Replace it if necessary
Dirt trap	Visual inspection	Every one month	Check and clean the foreign matter. Dry it by compressed air.

Table 3

8.2 Maintenance

8.2.1 Checking oil level

- 1) The oil level shall be always kept between position MAX and MIN of oil sight glass during operation (refer to Fig. 6). Add oil when the oil level is lower than position MIN and discharge oil by removing the oil drain plug (refer to Fig. 1) when the oil level is above position MAX.
- 2) Check the oil color. Replace the oil if the color is other than clear and transparent.

8.2.2 Checking pump's noise

The pump noise shall be continuous and stable during operation. If any abnormal noise is found, refer to Table 4 Troubleshooting.

8.2.3 Changing oil

- 1) Change the oil when the pump is switched off and cooled down.
- 2) Remove the oil drain plug (Fig. 1) and drain the used oil into a suitable receptacle. When the flow of oil stops, screw on the oil drain plug and start the pump shortly (10 seconds). Switch off the pump, remove the plug again and drain the residual oil. Screw on the oil drain plug (Check O-ring, replace it if necessary). Remove the oil fill plug (Fig. 1) and fill in fresh pump oil which is specified or provided by the manufacturer.
- 3) After oil changing, handle the used oil stored in an assigned container according to local relevant environmental laws and regulations.

8.2.4 Regularly checking intake filter and anti-suckback valve

Regularly take down the intake port and clean the filter by compressed air. Check the seal position of anti-suckback valve for cleanness, damage, hardening and height.

8.2.2 Checking pump's noise

Regularly remove the clean the fan cover and motor fan. Clean them by compressed air and reinstall them property before pump operation.

9. Troubleshooting

Fault	Possible reason	Remedy
Pump cannot start	Wiring is malfunctioning.	Check and repair wiring.
	Operation voltage is abnormal.	Make sure the voltage is within rated voltage's tolerance of $\pm 10\%$.
	Motor is malfunctioning.	Replace the motor.
	Overload protector is active.	Check the ambient temperature.
	Oil temperature is below 10°C .	Increase the ambient temperature.
	Pump is jammed.	Repair the pump.
	Out of operating for long.	Repair the pump.
	Oil is too viscous.	Change the oil.
	Exhaust filter or exhaust line is clogged.	Replace the filter or clean the exhaust line.
	Pump inner parts are damaged.	Repair the pump.
Pump cannot reach ultimate pressure.	Vacuum system configuration is unreasonable.	Choose a suitable pump.
	Vacuum system leak.	Check the system.
	Measuring technique or gauge is unsuitable.	Use correct measuring technique and gauge. Measure the pressure directly at pump's intake port.
	Oil filter is obstructed.	Change the oil filter.
	Oil is unsuitable.	Change the oil.
	Oil channel is obstructed.	Clean the channel.
	Oil is insufficient.	Add the oil to the level as requested.
	Anti-suck back oil valve is malfunctioning.	Repair anti-suck back oil valve.
	Intake line is dirty.	Clean vacuum lines.

Fault	Possible reason	Remedy
Pumping speed is too low.	Intake port line is clogged.	Clean the intake port line.
	Connecting lines are too narrow or too long.	Use adequately wide and short connecting lines.
	Exhaust line is clogged.	Keep exhaust port line unobstructed.
	Oil mist filter is clogged.	Change the oil mist filter.
	Anti-suck back oil valve is malfunctioning.	Repair anti-suck back oil valve.
Noise is abnormal	Operation voltage is abnormal.	Check the power supply, switches and wiring connection. Make sure the voltage is within rated voltage's tolerance of $\pm 10\%$
	Foreign matters entering into the pump.	Repair the pump.
	The oil level of the pump is too low.	Add the oil to the level as requested.
	The pump inner parts are damaged.	Disassembly and replace the parts.
Pump gets hotter than usually observed.	Poor ventilation.	Improve ventilation environment.
	The fan is damaged.	Change the fan.
	Pumped gas temperature is too high.	Add cold trap at intake port.
	The oil filter or exhaust line is clogged.	Replace oil filter or clean the exhaust line.
	Oil is unsuitable.	Change the oil.
	Oil channel is obstructed.	Clean the channel.
	Oil is insufficient.	Add the oil to the level as requested.
	The condenser is dirty.	Clean the condenser.
	Ambient temperature is too high.	Reduce ambient temperature.

7. Warranty

SH-VDC series vacuum pump have one year guarantee from the date of purchase. Spare parts will be provided free of charge within the period of guarantee in condition that the pump is operated according to the operating manual.

The following failures will be charged for repair:

- 1) Malfunction caused by nature disasters or artificial factors.
- 2) Malfunction caused by special operating environment.
- 3) Damage of wear parts.
- 4) Malfunction caused by abnormal operation or incorrect use indentified by our engineers.



Head office

www.silhumsil.com

Sejong, Republic of Korea

Tel. +82 44 866 5172 Fax. +82 44 866 5174

E-mail: sales@samheung21.com

Overseas sales office

www.shscientific.com

Daejeon, Republic of Korea

Tel. +82 42 826 5172 Fax. +82 42 826 5173

E-mail: i-sales@samheung21.com

SH Scientific Corporation USA

www.labandfurnace.com

Portland, Oregon, USA

E-mail: jhbang@labandfurnace.com